

Cat. No.
TLP-T01

ProNA™ Tension-Gel Making Solution

■ TLP-T01 0.75ml

This product is for laboratory research only and not for diagnostic use

Rn 202, 336-4 Jangdae-Dong, Yuseong-gu, Daejeon, Korea
Tel. 042-825-6992 Fax. 042-367-6991
E-mail: translab001@hanmail.net
Web: www.translabbio.co.kr

TransLab
Bioscience in KOREA

Description

- ProNA™ Tension-Gel Making solution can easily make Tension Acrylamide Gel if this is added to 30% acrylamide/bis solution.

Content

- Tension-Gel Making solution 0.75 ml
- User manual 1 Sheet

Applications

- Western blotting

Storage Information

- Store at RT

Protocol

- Tension-Gel Making Solution is added to 30% Acrylamide/Bis stock solution in a final concentration of 0.15%

0% Acrylamide/Bis Stock Solution	Tension-Gel Making Solution
500 ml	0.75 ml
250 ml	0.375 ml
100 ml	0.15 ml

Troubleshooting

- Please E-mail us at translab001@daum.net to request full detailed questions

Related Products

- 10x Effect running buffer	1 L	Cat. No. TLP-105
- 10x Cooling-Free transfer buffer	1 L	Cat. No. TLP-110
- Prestained protein marker	500 ul	Cat. No. TLP-111
- ECL Ottimo kit	500 ml	Cat. No. TLP-112.1
- Fast staining solution	500 ml	Cat. No. TLP-114
- 5X Phospho-block	500 ml	Cat. No. TLP-115.1P
- 5X General-block	500 ml	Cat. No. TLP-115.1G
- Western blot reprobe solution	500 ml	Cat. No. TLP-116
- CET1 lysis buffer	100 ml	Cat. No. TLP-121CETi